



How to use nonstandard filters

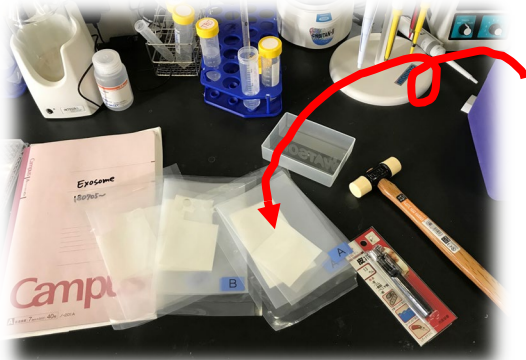

In our co-culture vessel, a circular filter with a diameter of 13 mm is standard, but if you want to use a filter other than 13 mm, adjust the size using a punch. Please purchase a commercially available punch 13 mm diameter. Please use Punch after cleaning and alcohol sterilization.

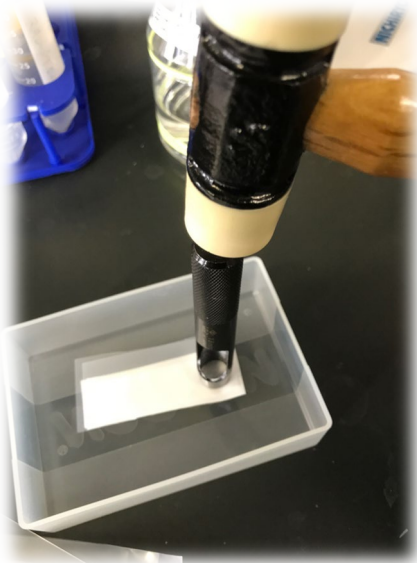

Take care not to scratch the experiment desk even if it punches out too much using the lid of the chip holder etc. Punch out with a hammer. First you need a knack, but it is easy if you get used to it.



After punch, the filter will be the perfect size. Please sterilize the filter with 70% ethanol.

